



## Indicator Status

Color	Status	Meaning
Green	Normally ON	Laser preheat indication, indicating that the laser is about to launch.
	Flash	—
Blue	Normally ON	Laser is emitting.
	Flash	Warning: Laser internal temperature exceeds 55 degrees Celsius. The laser remains emitting. Note: The flash will stop automatically when temperature will drop to 55 degrees Celsius.
Orange	Normally ON	Circuit fault, laser stops emitting.
	Flash	Warning: Laser internal temperature exceeds 85 degrees Celsius. The laser remains emitting. Note: The flash will stop automatically when temperature will drop to 55 degrees Celsius.
Red	Normally ON	Input voltage range overvoltage or undervoltage, laser stops emitting.
	Flash	—