


# Operation Instruction of PGL-III portable Laser

## 1. Opening method

- (1).Unscrew the end cap and install two “C” batteries
- (2).Screw the shutter of the laser.
- (3).Insert remote controller.

Note: Pay attention to remote controller notch. Insert tightly. Avoid damaging the laser.

- (4).Insert the key switch, push and screw the key switch about 90 degree clockwise.
- (5).Push switch button.

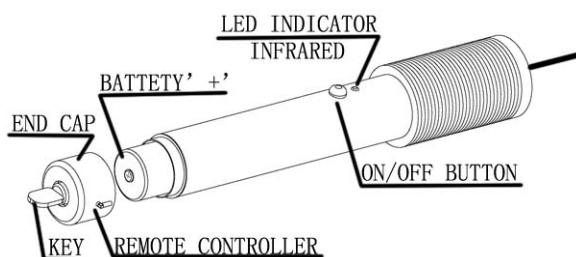

## 2. Closing method is reverse opening method.

## 3. The product is warranted for half year.

## 4. Attention: CNI does not assume liability for its laser products if any one of the following circumstances occurs.

- (1) Disassembly of the laser head by the user.
- (2) Influence to the proper laser performance as a results of injury to the outward surface.
- (3) No attachment of the original marks, e.g., serial numbers, of the laser head.
- (4) Invalidity of the repair guarantee after the specified warranty period.

## 5. Note

- (1).Laser is harmful to human being, and so you must avoid to be radiated directly. Especially, the laser beam should never point to eyes.
- (2).The laser should be operated in dry and clean environment.
- (3).Forbid shocking and oscillating.